

Manganese Field Test Kit

(Read Instructions before use)



ASIANMEDIC CO., LTD.

Tel. 6689-185-8999, 8890-898-5188

E-mail: sales@asianmedic.com



Procedure

Step 1

Add water sample into the reaction bottle to the level of the black line.



Step 2

Add reagent **Mn1** into the reaction bottle to the level of the red line. Swirl to mix.



Step 3

Add 2 spoons of reagent **Mn2** into the reaction bottle.





Step 4




Swirl to dissolve and wait for 10 minutes.

Step 5

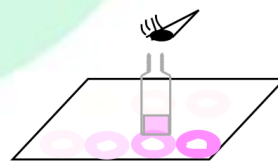Uncap and compare the developed color with the standard color scale by placing the uncapped reaction bottle at the center of color ring to match the color of the solution and the ring

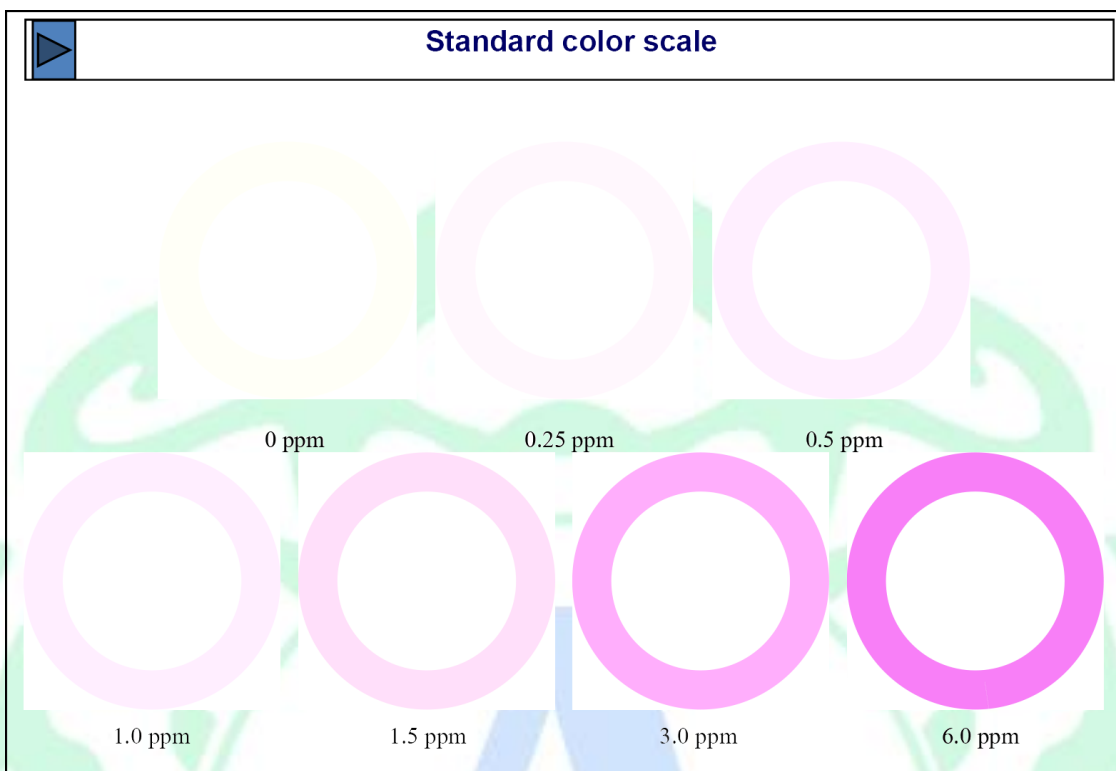



 Package contents	 Cautions
<ol style="list-style-type: none">1) Reaction bottle2) Syringe 5 ml3) Reagent Mn14) Reagent Mn25) Spoon (for handling reagent Mn2)	<ol style="list-style-type: none">1) Immediately close all reagent containers after use.2) Avoid contact of liquid reagents with skin and eyes.3) Always wash your hands after using the kit.4) Keep the kit out of direct sunlight and away from heat, food, pets and the reach of children
 Information on chemicals	
<p><u>Reagent Mn1</u> contains corrosive reagent that can cause severe burns to skin and eyes. Harmful if swallowed</p>	