

Nitrite Field Test Kit

Read instructions before use



ASIANMEDIC CO., LTD.

Tel. 6689-185-8999, 8890-898-5188

E-mail: sales@asianmedic.com



Package contents

- | | |
|------------------------------|----------|
| 1. Reaction bottle | 2 piece |
| 2. Syringe 1ml | 1 piece |
| 3. Reagent N 1 | 1 bottle |
| 4. Reagent N 2 | 1 bottle |
| 5. Plastic bag (paper strip) | 5 piece |



Cautions

- 1) Immediately close all reagent containers after use.
- 2) Avoid contact of liquid reagents with skin and eyes.
- 3) Always wash your hands after using the kit.
- 4) Keep the kit out of direct sunlight and away from heat, food, pets, and the reach of children.



Information on chemicals

Reagent N1 contains corrosive reagent that can cause severe burns to skin and eyes. Harmful if swallowed.



Testing Procedure

Step 1

Add 1 ml of water sample into the reaction bottle1 with syringe 1ml-contained-syringe



Step 2

Add 1 drop of reagent N1 and swirl gently to mix the solution



Step 3

Immerse the strip N2 into the solution



Step 4

Compare the developed colour with the standard colour scale

สีที่ปรากฏ								
ความเข้มข้นของไนไตรท์ (ppm N)	0.0	0.01	0.05	0.1	0.2	0.5	1.0	2.0



Standard color scale

Color Scale								
Concentration of nitrite (ppm N)	0.0	0.01	0.05	0.1	0.2	0.5	1.0	2.0